

Additional Vial Instructions

Additional vials may be added to any ProEdge Dental Water Labs testing kits.

1. Label Vials

With a permanent marker, number each additional vial sequentially, starting with the next available

number. For example, if you ordered a 6-Vial Test Kit, label your additional vial 7 and then 8 and so forth.



2. Collect Samples

Fill sample vial to 90% capacity. Close lid securely.



FLIP FOR STEP 3



3. Complete Submission Form

Fill out the form **completely** to ensure accurate results. The vial number you put on the sample will be indicated on your results.



ProEdge
DENTAL WATER LABS

We'd love to help.

888.843.3343

ProEdgeDental.com